



MDS-15

High-throughput Microwave Sample Preparation Workstation

Fully Functional, Multiple Use, Safe Operation



The Transformer of Microwave Digestion

MDS-15

*High-thro
SO*

Fully Functional, Multiple Use, Safe Operation

The Transformer of Microwave Digestion

- The quartz digestion vessel enables the digestion process to be visible;
- Initiative wireless remote video monitoring system, together with quartz vessels, makes the reaction be seen clearly at a glance. Pictures and videos can be taken during the entire process;
- 500 ml ultra-large digestion vessel meets the special requirements for the digestion of samples of up to 10 grams;
- The drying rotor aides in the drying process of samples;
- Sineo's patented safety bolt design eliminates the need for explosion-proof membranes and other costly consumables;
- The aerospace composite fiber outer vessel ensures the highest level of vessel security;
- Large 65L stainless steel chamber and dual magnetron structure, is designed for high energy and uniform microwave field;
- With over 20 years of experience in the industry Sineo is known as an innovation leader with numerous patents;



High-Throughput Microwave Sample Preparation Workstation

Our Latest microwave system (The MDS-15) is the culmination of over 20 years of refinement. The MDS-15 is a high-throughput, multi-functional closed microwave sample preparation workstation with the latest technology and a lifetime warranty on the core component. **Versatility:** The MDS-15 is a versatile workstation which can handle the routine work, and more complicated and demanding samples in your prep rack with a wide variety of vessels to choose from. **Safety is priority number one!** At Sineo we put safety at the top priority, which is why we have painstakingly designed every aspect of our microwave system with the highest level of security measures, such as the use of aerospace composite fiber vessel sleeves, and our patented Safety Bolt. **Efficient by design:** MDS-15 applies high-energy uniform microwave field, it delivers a rapid digestion with a quick 20 minute cool down time. Sineo has incorporated the latest technology of material design while engineering a very safe, durable and easy to use microwave.



Functional vessels for your various application needs

Conventional digestion

With 16 high-throughput GP-100 digestion vessel set, MDS-15 can work for long time digestion at ultra high temperature. It meets the requirements for most of the digestion tasks; the new 8 QZ-100 quartz closed vessel set dramatically improves the digestion capability and visibility, even the most demanding materials could be successfully digested. With remote Wi-Fi video monitoring system for quartz vessel, the whole process can be observed, pictures and videos are available during the entire process.

Ultra-high sample amounts digestion

LV-500 digestion vessel carousel enables eight 500ml ultra large vessels to digest simultaneously, upgrading the sample amount to an unparalleled level of 10 gram. With magnetic stirring feature, sample digestion could be done more quickly and thoroughly. Innovative Safety Bolt design ensures the combination of vertical & quantitative pressure relief and complete digestion. The digestion tasks of high sample amount thus could be accomplished.


Sample drying

The F-DRY drying rotor supports rapid and uniform microwave drying, the sample drying time is only a quarter that of traditional method, it can also prevent sample contamination and carbonation. There' s no need of any modification to the main machine to complete the drying task.

Microwave extraction, microwave Kjeldahl Nitrogen Determination, etc

Sineo' s application engineers have developed the application methods for microwave extraction, microwave Kjeldahl nitrogen determination, etc. Our technical support is at your service all the time.




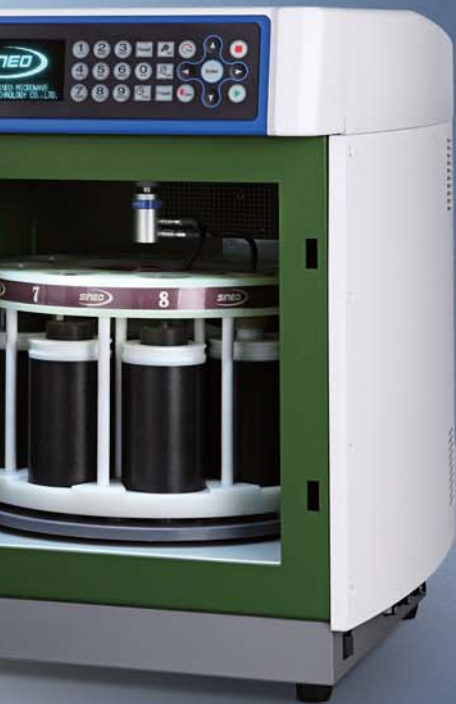

- The outer vessel of MDS-15 exclusively made by ultra-strength aerospace composite fiber is invincible in anti-explosion, and its performance indicators such as corrosion resistance, high temperature/impact/pressure resistance are far better than that of the widely used modified PEEK engineering plastics vessel (this material is fusible at high temperature, fragile at high pressure and explosive by chemical corrosion), fundamentally eliminating safety risks to operator in use.



- Quantified safety bolt design, ensures samples be closed completely and triggers a quantified pressure relief while over pressure; Safety Bolt unit, instead of safety membrane and other consumables, ensures the digestion vessel be sealed completely under normal working conditions, which avoids the loss of volatile elements and incomplete digestion. And only when the pressure is too high and may constitute a danger to the safety, the safety bolt will automatically blow out vertically and the cover auto-up to release the pressure, achieving quantified vertical blast pressure-relief to guarantee its well operation. Under normal operation, the safety bolt requires no replacement. In addition, it is easy for venting to open the cover after completion of digestion.

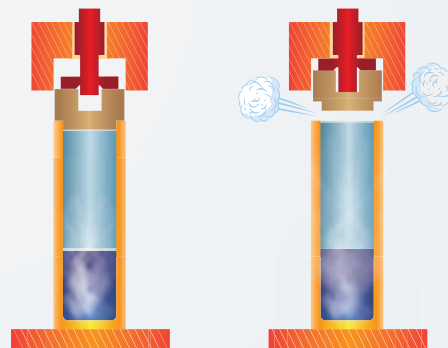

- The industrial-leading pressure measuring technology by piezoelectric crystal and high-precision Pt sensor temperature measurement, through closed-loop control of microwave power by inverter technique, ensure the accuracy of pressure and temperature monitoring and control. The application of patented piezoelectric crystal brings complete isolation of samples from pressure measurement system in digestion process, thoroughly solving the problems of cross contamination of samples due to commonly used air pipe in market and of the limitation in digestion samples because of low-pressure proof of air pipe.



- The whole set of digestion vessels was designed to rotate continuously in one direction, it breaks conventions of $<360^\circ$ back and forth rotation, avoiding uneven heating on vessels and reducing impact on turntable motor to extend service life. The temperature and pressure cables are short, that they are free from twisting in running and can be disassembled easily.
- The new industrial design is smart and colorful, that in line with the needs of modern laboratory construction needs. It re-layout its internal structure more scientifically, with the 65L industrial 316L stainless steel reactor chamber and dual magnetron structure, ensures uniform heating of the microwave and high microwave power emitting.





- Innovative remote wireless video monitoring system, together with the high-intensity transparent quartz vessel, makes the digestion process clearly visible. Mobile phones and tablets can realize the observation and monitoring of the reaction, making the operation safer and more convenient.



- Sturdy and durable industrial-grade chamber design strengthen its impact resistance, and multi-layer chemical resistant coating greatly improves the service life and safety of the system; the popup cushioning explosion-proof sliding chamber door (Auto-Pop structure) builds a passive safety protection system, and it can disable the microwave power automatically once explosion happens; efficient exhaust system design achieves fast and safe forced air cooling (20 min cooling time from 200°C to 60°C), improving operational efficiency.



The Popup Cushioning Explosion Proof Sliding Chamber Door

High-performance digestion rotors



16 GP-100 ultra-high-strength frame closed vessel

16 high-throughput GP-100 digestion vessel set adopts the latest technologies and high-end materials, it is capable of long time digestion work under high temperature and pressure, which can meet the needs of vast majority of digestion tasks. It is especially suitable for the test house or labs handling with various sample types and large capacity. GP-100 is also suitable for the applications of solvent extraction, microwave kjeldahl nitrogen determination and microwave hydro-thermal synthesis, etc.

| | |
|------------------------------|--|
| Maximum design pressure: | 10MPa (1500psi) |
| Maximum design temperature: | 300°C |
| Internal vessel volume: | 100ml |
| Outer vessel material: | Ultrastrength aerospace composite fiber, high temperature and pressure resistance reach 600°C and 66 Mpa (10000psi). |
| Inner vessel material: | TFM (modified polytetrafluoroethylene) |
| Maximum processing quantity: | 16 vessels |

8 QZ-100 high temperature/pressure visual closed quartz vessel

8 QZ-100 quartz digestion vessel set is developed by Sineo according to market demand and the latest technology in the industry. It is able to work at even tougher condition than GP-100, under higher temperature and pressure, even the most demanding material can be successfully digested. At the same time, the remote video monitoring system provided can carry out real-time monitor of the reaction process inside the vessel. Mobile phone or tablet with monitoring App will help the operator to follow up the digestion reaction and enjoy a safer operation.

| | |
|------------------------------|----------------------------|
| Maximum design pressure: | 10MPa (1500psi) |
| Maximum design temperature: | 310°C |
| Internal vessel volume: | 100ml |
| Inner vessel material: | ultra-pure quartz material |
| Maximum processing quantity: | 8 vessels |
| Resistance to HF: | No |





8 LV-500 ultra large capacity digestion vessel carousel

LV-500 digestion vessel carousel provides 8 ultra-large vessels with 500ml in volume, the sample amount can be increased to unparalleled 10 gram grade. For testing of trace element, there is no need to collect samples for many times. It improves efficiency and at the same time avoids the operation error, thus the needs for high sample amount digestion can be achieved. At the same time, the full vessel magnetic stirring helps the sample to digest quickly and thoroughly. The aerospace fiber outer vessel, high-strength vessel frame and Safety Bolt design, ensure the safe, vertical and quantitative pressure relief as well as the thorough digestion.

| | |
|------------------------------|---|
| Maximum design pressure: | 4Mpa (600psi) |
| Maximum design temperature: | 230°C |
| Internal vessel volume: | 500ml |
| Outer vessel material: | Ultrastrength aerospace composite fiber, high temperature and pressure resistance reach 600°C and 66 Mpa(10000psi). |
| Inner vessel material: | TFM (modified polytetrafluoroethylene) |
| Maximum processing quantity: | 8 vessels |

F-DRY microwave drying rotor

F-DRY microwave drying rotor realizes rapid and uniform drying, it takes the moisture off sample through microwave heating and vortex air. The sample drying time is only a quarter of that required by conventional heating. The special design of the air inlet and outlet can prevent sample from contamination and loss, at the same time, there is no carbonized risks of the sample by microwave heating. The heating task can be completed without any modification of the main machine. It meets the standard for special sample testing like soil.






| | |
|------------------------|---|
| Size of drying plate: | 27cm in diameter |
| Drying plate material: | TFM (modified polytetrafluoroethylene) |
| Drying wind speed: | three adjustable grades of high, middle and low drying wind |

MDS-15 High-throughput Microwave Sample Preparation Workstation

Main Technical Parameters:

| | |
|--|---|
| Power | 220~240 VAC 50/60Hz 16A |
| Microwave source | 2450MHz, dual magnetron design |
| Installed power | 3600W |
| Maximum output power | 2200W, non-pulse continuous automatic variable frequency control |
| Cavity | 65L industrial grade 316L stainless steel chamber, multi-layer chemical resistant PFA coating |
| Door design | 6 layer steel structure and popup cushioning explosion-proof sliding chamber door (Auto-Pop) design, electronic and mechanical dual-control door lock |
| Video monitoring system (Optional) | Wireless video monitoring system, mobile phones and tablets can realize the observation and monitoring of the reaction |
| Pressure measurement and control system | Piezoelectric crystal pressure sensor, pressure control range :0~10MPa (1500 psi), accuracy ± 0.01 MPa |
| Temperature measurement and control system | High-precision platinum resistor temperature sensor, temperature range :0-300°C, accuracy $\pm 1^\circ\text{C}$; IR temperature sensor as optional, temperature range :0-300°C, accuracy $\pm 1^\circ\text{C}$. |
| Outer vessel material | Explosion-proof outer vessel made of aerospace composite fiber. high temperature and pressure resistance reach 600°C and 66 Mpa(10000psi) |
| Inner vessel material | Modified TFM material |
| Chamber exhaust system | High-power anticorrosion axial fan,20 minutes cooling time (from 200°C to 60°C) |
| Working environment | 0~40°C, 15~80%RH |
| Whole physical size/weight | 550 × 670 × 700mm, 70 KG |

| Name of reaction vessel | 16 GP-100 ultrahigh strength frame closed vessel | 8 QZ-100 high temperature/pressure visual closed quartz vessel | 8 LV-500 ultra large capacity digestion vessel carousel |
|---------------------------------------|---|--|---|
| Model of reaction vessel | GP-100 | QZ-100 | LV-500 |
| Processing capacity in one batch | 16 | 8 | 8 |
| Inner vessel material | TFM | Pure Quartz | TFM |
| Explosion-proof outer vessel material | Aerospace composite fiber Xtra Fiber | N/A | Aerospace composite fiber Xtra Fiber |
| Frame structure of digestion rotor | Independent frame structure of each digestion vessel | Independent frame structure of each digestion vessel | Integrated high strength alloy rotary carousel |
| Reaction vessel volume | 100ml | 100ml | 500ml |
| Maximum design pressure | 10MPa(1500psi) | 10MPa (1500psi) | 4MPa (600psi) |
| Maximum temperature | 300°C | 310°C | 240°C |
| Maximum sample amount | 0.5-1.0g | 0.5-1.0g | 5-10g |
| Resistance to HF | Yes | No | Yes |
| Magnetic stirring function | No | No | Yes |
| Typical application | Food, soil, metal, alloy, cosmetic, plant, etc. | Fat, oil, plastic, coal, etc. | Biological samples, environmental samples, etc. |
| Photo |  |  |  |

Fully Functional, Multiple Use, Safe Operation

The Transformer of Microwave Digestion

